

O-Mount

(An organic based mounting media for immunohistochemistry)

Storage: Room Temp.
Do Not Freeze.

Catalog No.: E02-18 18ml

Intended Use:

O-Mount is designed as a routine coverslipping mounting media for organic insoluble chromogens, such as DAB. It also works well for routine H & E staining. O-Mount is soluble in both xylene and toluene. O-Mount is supplied in a 18 ml glass dropper bottle.

Kit Components:

Product	Catalog No.	Liquid media
O-Mount	E02-18	18 ml

Recommended Protocol:

1. Remove slides from clearing reagent, shake off excess liquid from the slide and quickly wipe the back of the slide.
2. Invert dropper bottle to displace bubbles from the tip-end.
3. Squeeze out the first drop onto a paper towel to remove of air bubble on the tip.
4. Add one or two drops of O-Mount on the tissue section.
5. Carefully lay down coverglass to cover the tissue section
6. Gently press out any noticeable air bubbles
7. Use the minimal amount of media necessary to effectively bond the coverglass to the slide
8. Allow sufficient time of drying slides before storage.

Precautions:

O-mount is flammable. Store in containers or area that is fire prove. Do no store near heat, sparks, flames or strong oxidants. Keep bottle capped when not in use.

Remarks:

For research use only. Not for diagnostic or therapeutic use.